

H&E Staining Protocol

Rehydration

- Xylene 5min 2X
- 100% EtOH 1min
- 95% EtOH 1min
- 80% EtOH 1min
- 70% EtOH 1min
- dH₂O

Staining

- Hematoxylin 1-5min
- Rinse dH₂O
- Acid Alcohol 2sec. (.1% HCL in 50% EtOH)
- Rinse Tap Water 15min
- Eosin 1min
- Rinse in a few quick changes of dH₂O

Dehydration

- 95% EtOH about 1 min as needed
- 100% EtOH 1min
- Xylene 5min 2X
- Coverslip